

# Laboratory protocol: isolation of rat liver mitochondria

Sumbalova Z<sup>1,2</sup>, Fontana-Ayoub M<sup>2</sup>, Krumschnabel G<sup>2</sup>

<sup>1</sup>Pharmacobiochem Lab, Fac Medicine,  
Comenius Univ, Bratislava, Slovak Republic

<sup>2</sup>**Oroboros Instruments**

O2k High-Resolution FluoRespirometry (HRFR)  
Schöpfstr 18, A-6020 Innsbruck, Austria  
Email: [instruments@orooboros.at](mailto:instruments@orooboros.at)  
[www.orooboros.at](http://www.orooboros.at)

## 1. Preparation

Switch on centrifuge and let it cool down to 4 °C. Keep all buffers, dissection gear, homogenization tools and centrifuge rotor at 4 °C (or on ice).

### 1.1. Anesthesia

Rats are anesthetised by intraperitoneal injection of Thiopental (0.1 g/kg) or by CO<sub>2</sub> narcosis.

### 1.2. Isolation procedure

1. Kill rat, dissect out liver (take weight) and put liver in ice-cold isolation medium.
2. Determine wet weight, take 1.5 g of the tissue for isolation.
3. Transfer the tissue to a pre-cooled glass beaker (20 ml) with ice-cold isolation medium, discard all medium.
4. Mince the tissue into small pieces using a pair of sharp scissors (should become a mash), add drops of medium while cutting.
5. Suspend with ~ 5 - 10 volumes of ice-cold isolation medium and transfer to a pre-cooled glass/Teflon potter.
6. Homogenize the tissue with 8 - 10 strokes at 1,000 rpm, add more media.
7. Transfer to the 50 ml Falcon tube, bring the volume to get <10% homogenate (1 g tissue to 15 - 20 ml homogenate).
8. Centrifuge at 1,000 *g* for 10 min at 4 °C.
9. Filter the supernatant through two layers of gauze (optional), transfer it into a new tube and centrifuge at 6,200 *g* for 10 min at 4 °C.
10. Discard the supernatant and re-suspend the mitochondria (sediment) in 15-20 ml of isolation medium per g tissue, centrifuge at 6,200 *g* for 10 min at 4 °C.
11. Discard the supernatant and re-suspend mitochondria in a small volume of the medium (the volume of mitochondrial suspension from 1.5 g tissue ~ 1.5 ml).
12. Store mitochondria on ice, use within 3-4 h.
13. Transfer subsamples (20 µl) into Eppendorf tubes and store at -20 °C for further analysis (protein concentration, citrate synthase).

## 2. Media

### 2.1. Isolation buffer

Chemical	Final conc. [mM]	Add to 1000 ml [g]
Mannitol	225.0	40.99
Succrose	75.0	25.67
EDTA	0.2	0.0744

Adjust pH to 7.4 with Tris, HCl.

## 3. References

Isolation protocol modified after Kucharska et al 2000:

Kucharská J, Braunová Z, Ulicná O, Zlatos L, Gvozdjáková A (2000) Deficit of coenzyme Q in heart and liver mitochondria of rats with streptozotocin-induced diabetes. *Physiol Res* 49:411-8. »[Bioblast link](#)«




O2k-Protocols

[http://wiki.orooboros.at/index.php/O2k-mitochondrial\\_preparations](http://wiki.orooboros.at/index.php/O2k-mitochondrial_preparations)

## Acknowledgements

Sumbalova Z was supported by Action Austria – Slovakia (2015-Feb-01 to Jun 30). Protocol development supported in part by K-Regio project **MitoFit**.  
[www.mitofit.org](http://www.mitofit.org)

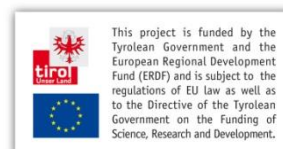

This project is funded by the Tyrolean Government and the European Regional Development Fund (ERDF) and is subject to the regulations of EU law as well as to the Directive of the Tyrolean Government on the Funding of Science, Research and Development.